

## UL-SB0012 New Control Computer Setup

### Tool List

1. New control computer with 64-bit Windows 10.

### Instructions

1. NOTE: Do not connect the computer to the UltraLo-1800 until instructed to do so as this may cause driver errors.
2. Complete the Windows installation process as per your organization's typical standards.
  - a. It's highly recommended that the user account that controls the UltraLo-1800 be named "UltraLo", and be the primary (preferably sole) account on the computer. Doing otherwise may cause errors.
  - b. It's recommended (but not required) that the computer be named "UL-[4-digit serial number]", eg "UL-1234".
  - c. If you wish to use the remaining standard settings XIA uses, please contact us at [support@xia.com](mailto:support@xia.com), but note that these settings change frequently with Windows versions so XIA's instructions may be out of date.
3. Disable automatic restart for Windows Updates via group policy.
  - a. Open the start menu, search for "group policy".
  - b. Navigate to Local Computer Policy -> Computer Configuration -> Administrative Templates -> Windows Components -> Windows Update.
  - c. Double-click on "Configure Automatic Updates"
    - i. Select the "Enable" radio button.
    - ii. Set "Configure automatic updating" to "3"
    - iii. Press "OK"
  - d. Double-click on "No auto-restart with logged on users for scheduled Automatic Update installation"
    - i. Select the "Enable" radio button.
    - ii. Press "OK".
  - e. Close the group policy window.
4. Disable sleep/hibernation
  - a. Open the Control Panel (Start-> search for "Control Panel").
  - b. Navigate to Power Options and click "Change plan settings" for the selected plan.
  - c. Set the "Sleep" option to "Never" (this may require administrator permissions).
  - d. Click "Change advanced power settings".
  - e. Expand all the menus under "sleep" (eg "Sleep After", "Hibernate After") and ensure they're all turned off.
5. Download the latest CounterMeasure release from <http://support.xia.com/default.asp?W430> and install it.

6. Download the latest version of the SeaMAX program for the optical isolator from <https://www.sealevel.com/support/software-seamax-windows/> and install it.
  - a. Note: installation may prompt for the product name, it's "8112".
7. Download and install the serial power driver from <https://www.advantech.com/support/details/driver?id=1-13U9QTV>.
8. Connect the computer to the UltraLo-1800 and turn on the power.
9. Start CounterMeasure.
  - a. CounterMeasure doesn't start, contact XIA at [support@xia.com](mailto:support@xia.com) and include the "CounterMeasure.log" file found in C:\Users\UltraLo\AppData\Roaming\XIA LLC\CounterMeasure\Logs, as well as a picture of the device manager with the "Universal Serial Bus controllers" and "Ports (COM & LPT)" sections expanded.
10. Update the calibration certificate with the one provided by XIA.
  - a. Restart CounterMeasure for all the settings in the certificate to take effect.